
TOPSPIN Control and Function Keys

Focus anywhere in TOPSPIN

Esc	Put the focus in the command line
Shift+Esc	Display menu bar and toolbars (if hidden)
F2	Put the focus in the browser / portfolio
F1	Search string in command help or NMR Guide [help]
F6	Select the next window in the data area
Alt+F4	Terminate TOPSPIN [exit]
Ctrl+d	Switch the browser/portfolio on/off
Ctrl+o	Open data [open]
Ctrl+f	Find data [find]
Ctrl+n	New data [new]
Ctrl+p	Print current data [print]
Ctrl+s	Save current data [sav]
Ctrl+w	Close active window [close]
Ctrl+c	Copy a text that you selected/highlighted in an error box, dialog box, pulse program, title etc., to the clipboard
Ctrl+v	Paste text from the clipboard into any editable field.

Focus in the Command Line

Ctrl+Backspace	Kill current input
Ctrl+Delete	Kill current input
UpArrow	Select previous command (if available).
DownArrow	Select next command (if available).

Focus in the Browser/Portfolio

UpArrow	Select previous dataset
DownArrow	Select next dataset
Enter	Browser: expand selected node
Enter	Portfolio: display selected data

Focus anywhere in TOPSPIN

<i>Scaling Data</i>	
Alt+PageUp	Scale up the data by a factor of 2 [*2]
Alt+PageDown	Scale down the data by a factor of 2 [/2]
Ctrl+Alt+PageUp	Scale up by a factor of 2, in all data windows
Ctrl+Alt+PageDown	Scale down by a factor of 2, in all data windows
Alt+Enter	Perform a vertical reset
Ctrl+Alt+Enter	Perform a vertical reset in all data windows
<i>Zooming data</i>	
Alt+Plus	Zoom in [.zi]
Alt+Minus	Zoom out [.zo]
Ctrl+Alt+Plus	Zoom in, in all data windows
Ctrl+Alt+Minus	Zoom out, in all data windows
<i>Shifting Data</i>	
Alt+UpArrow	Shift spectrum up [.su]
Alt+DownArrow	Shift spectrum down [.sd]
Alt+LeftArrow	Shift spectrum to the left [.sl]
Alt+RightArrow	Shift spectrum to the right [.sr]
Ctrl+Alt+UpArrow	Shift spectrum up, in all data windows
Ctrl+Alt+DownArrow	Shift spectrum down, in all data windows
Ctrl+Alt+LeftArrow	Shift spectrum to the left, in all data windows
Ctrl+Alt+RightArrow	Shift spectrum to the right, in all data windows

Focus in a Table (e.g. peaks, integrals, nuclei, solvents)

delete	Delete the selected entries
home	Select the first entry
end	Select the last entry
Shift+Home	Select the current and first entry and all in between
Shift+End	Select the current and last entry and all in between
DownArrow	Select next entry
UpArrow	Select previous entry
Ctrl+a	Select all entries
Ctrl+c	Copy the selected entries to the clipboard
Ctrl+z	Undo last action
Ctrl+y	Redo last undo action

Note that function keys can be redefined as described in the TOPSPIN Users Guide.

Focus in a Plot Editor

F1	Open the Plot Editor Manual
F5	Refresh
Alt+F4	Close the Plot Editor window
Ctrl+F6	Display next layout
Ctrl+Shift+F6	Display previous layout
Ctrl+tab	Display next layout
delete	Delete the selected objects
Ctrl+a	Select all objects
Ctrl+c	Copy the selected object from the Clipboard
Ctrl+g	Group the selected objects
Ctrl+Shift+g	Ungroup the selected object
Ctrl+i	Open TOPSPIN Interface
Ctrl+l	Lower the selected object
Ctrl+m	Unselect all objects
Ctrl+n	Open a new layout
Ctrl+o	Open an existing layout
Ctrl+p	Print the current layout
Ctrl+q	Close the Plot Editor window (Linux only)
Ctrl+r	Raise the selected object
Ctrl+s	Save the current layout
Ctrl+t	Reset X and Y scaling of all marked objects
Ctrl+v	Paste the object from the Clipboard
Ctrl+w	Open the attributes dialog window.
Ctrl+x	Cut the selected object and place it on the Clipboard
Ctrl+z	Undo the last action

